



# Veratox<sup>®</sup> for Hazelnut Allergen

**Note: Please read kit instructions completely before performing test.**

**Questions? Call 800/234-5333 or 517/372-9200**



1. Add 150  $\mu$ L controls and extracted samples to transfer wells.



2. Transfer 100  $\mu$ L to the antibody wells. Incubate for 10 minutes. Mix for 20 seconds by sliding back and forth on a flat surface.



3. Dump liquid from antibody wells.



4. Wash wells thoroughly. Repeat wash step 5 times with washing solution.



5. Tap out water on absorbent paper towel.



6. Transfer 100  $\mu$ L conjugate from reagent boat to antibody wells using 12-channel pipettor. Incubate for 10 minutes. Mix for 20 seconds by sliding back and forth on a flat surface.



7. Repeat steps 3–5 by dumping out the liquid, thoroughly washing the wells and tapping dry.



8. Transfer 100  $\mu$ L substrate from reagent boat to antibody wells using 12-channel pipettor. Incubate for 10 minutes. Mix for 20 seconds by sliding back and forth on a flat surface.



9. Transfer 100  $\mu$ L Red Stop from reagent boat to antibody wells. Mix.



10. Read results in a microwell reader.

## Product Specifications

**Format:** Sandwich ELISA

**Range of quantitation:** 2.5–25 ppm hazelnut

**Controls provided:** 0, 2.5, 5, 10 and 25 ppm

**Testing time:** 30 minutes

## Ordering Information

8420 Veratox for Hazelnut Allergen

Call **800/234-5333** to order or visit us online at [www.neogen.com](http://www.neogen.com)



800-234-5333 (USA/Canada) or 517-372-9200  
[foodsafety@neogen.com](mailto:foodsafety@neogen.com) • [foodsafety.neogen.com](http://foodsafety.neogen.com)

